



16. Back out and lower the Straddle forks and remove the Roller Truck from the straddle.
17. Position the Straddle and the LOS Transport Cart in close proximity to each other to facilitate easy transferring of the LOS to the Straddle Breadboard 1.
18. Place the *Teflon Sheeting* on the Optical Breadboard
19. Place the *Teflon Step Blocks* onto the Teflon sheet.

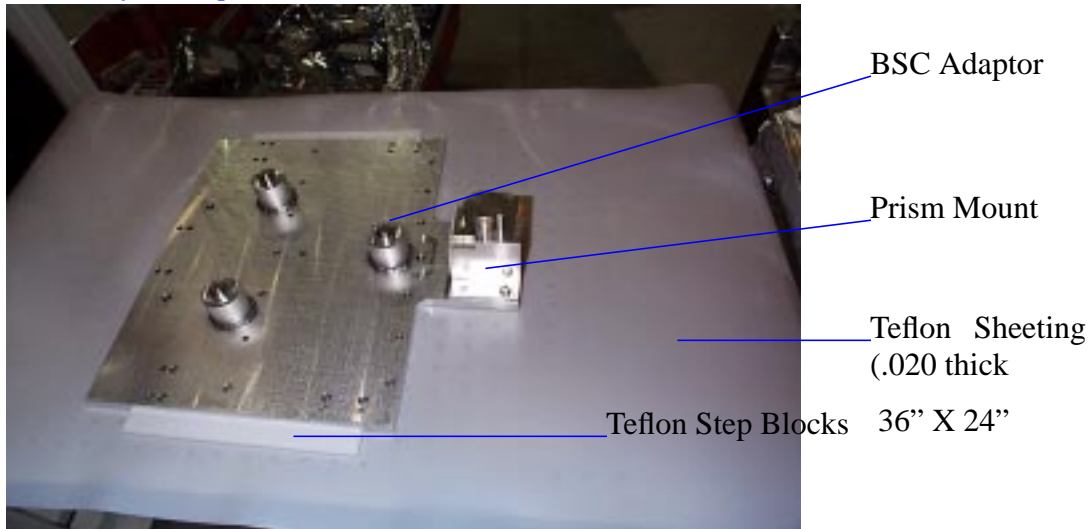

20. Place the *Alignment Fixture BSC Adaptor* onto the Teflon Step Blocks.

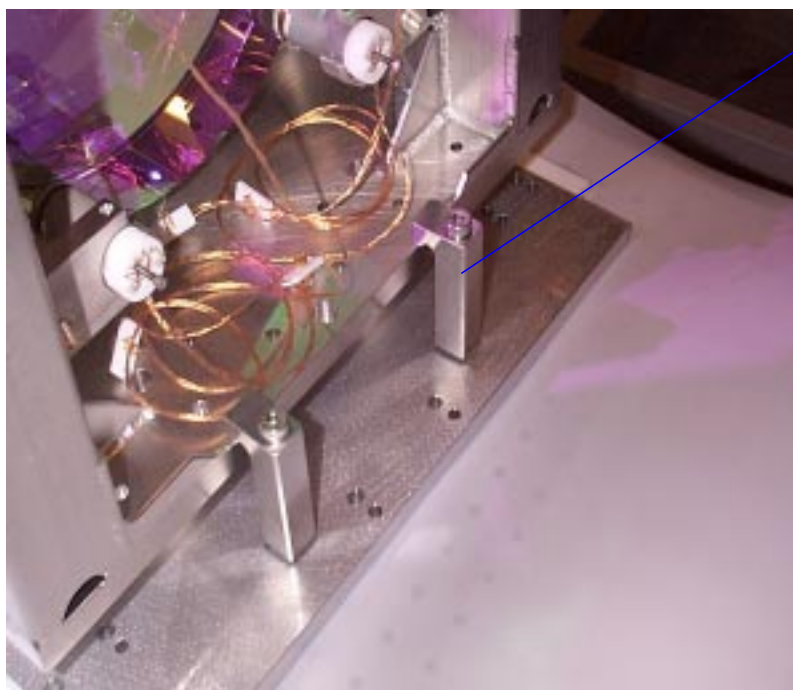

“Dog” clamps holding the structure to the BSC Adaptor Plate.

SUS cage with optic locate on the three posts and secured with the ‘dog’ clamps. Prism mount is on the front side (HR side of the optic)